# TOP NEWS

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***THE MOST ECO-FRIENDLY SAMPLE  
PREPARATION STATION!***

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Again and again media reports about the waste of resources and the increasing pollution with different gases. These are topics that concern all of us. We have taken these issues to heart and created a new solution that commits itself to **preserving energy and resources**.

Our latest sample preparation station is the **most eco-friendly special working table** by KUGEL medical for a variety of reasons!

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**MAKE A SUSTAINABLE CONTRIBUTION TO ENVIRONMENTAL PROTECTION WITH KUGEL MEDICAL**

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The **PEFC and FSC certificates** guarantee that the wood and raw materials we use for manufacturing our furniture comes from sustainably managed forests.

Additionally, for this sample preparation station we use an **eco-friendly permanganate carbon filter** which not only effectively vacuums away all harmful vapors and hence complies with the maximum concentration values at the workplace but is also economical in use and consequently **reduces your overheads by up to 30%**.

# THE MOST ECO-FRIENDLY SAMPLE PREPARATION STATION FROM KUGEL MEDICAL

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Our sample preparation station comprises of three parts that could, if necessary, also be obtained individually in the course of modernization works: **laboratory table with table plate made of high pressure laminate, perforated working area and active-carbon air recirculation system.**

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The laboratory plate is made of eco-friendly high pressure laminate. All raw materials and woods that are used during production come from **sustainably managed forests with PEFC- and FSC-certificates**. The surfaces can be individually color matched with any corporate identity. Additionally, the high pressure laminate plates are coated with the **antimicrobial Sanitized® coating with silver ions** which kills 99.9% of the bacteria within 24 hours. A C-shaped base frame is mounted under the table plate to provide plenty of leg room when working.

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**Moreover, high pressure laminate has a lot of outstanding features:**

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The perforated working area is inserted in the laboratory plate. All harmful vapors are being vacuumed downwards throughout the entire working area. When doing so, the **maximum concentration values for work involving formaldehyde are being underscored.**

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An active-carbon air recirculation system is placed under the laboratory table. The **integrated permanganate carbon filter** purifies the contaminated air and directs the clean air back to the laboratory. The great advantage in this connection is that a customer' supplied ventilation is not necessary. Thanks to its large permanganate carbon filter, the table can **be in use for up to 800\* working hours**. The filter can be easily replaced through the large inspection hole in the housing of the air recirculation system.

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\* operation time depends of various factors

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## Laboratory Table with Table Plate:

Material Table Plate: 38 mm high pressure laminate  
 Coating: antibacterial Sanitized®  
 coating with silver ions

## C-Shaped Base Frame:

Material: precision steel pipe as per  
 DIN 2395  
 Coating: powder coating  
 Color: light grey, RAL 7035

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## Perforated Working Area:

Material: stainless steel, material No. 1.4301  
 Perforation Working Plate: D17, T60  
 Air flow: min. 0.4 m/sec  
 Dimensions: 600 x 500 mm

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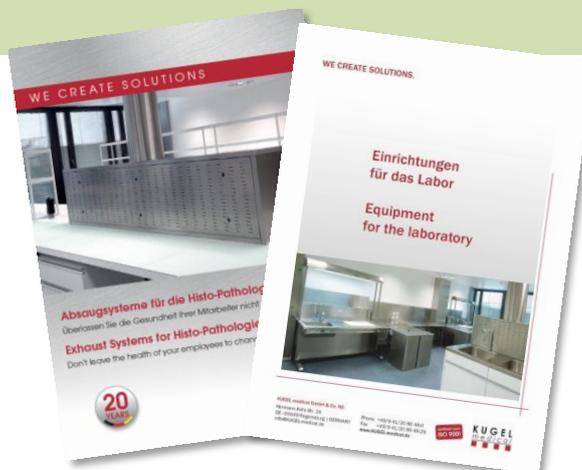
## Active-Carbon Air Recirculation System:

Material Housing: stainless steel  
 Working Temperature: max. +80 °C  
 Max. rel. Humidity: 70%  
 Max. Air capacity: approx. 350 m³/h  
 Delivery Includes: HEPA Filter H13,  
 Operating Hour Counter  
 Dimensions: 1000 x 385 x 455 mm

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**Individual solutions with customized dimensions  
 are available upon request!**

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**For more information, ask  
 now for our laboratory  
 furniture catalogue and  
 exhaust systems  
 brochure!**

**KUGEL medical GmbH & Co. KG**  
 Hermann-Köhl-Straße 2 A  
 DE – 93049 Regensburg  
 GERMANY

Phone +49 (0) 9 41/20 86 48-0  
 Fax +49 (0) 9 41/20 86 48-29  
 E-Mail info@KUGEL-medical.de  
[www.KUGEL-medical.de](http://www.KUGEL-medical.de)

**KUGEL**  
 medical